


# AppliChrom® ABOA SugarSep-Oligo-Ag

13062020 Application Note


## High Resolution HPLC Oligosaccharidanalytics in Water with RI-Detection

**AppliChrom ABOA SugarSep-Oligo-Ag, 300x8mm, Order No. SASOLAg103008**  
 "Maltodextrin 12 + extra glucose", Eluent: H<sub>2</sub>O, 0,25ml/min, 80°C



Total chromatogram

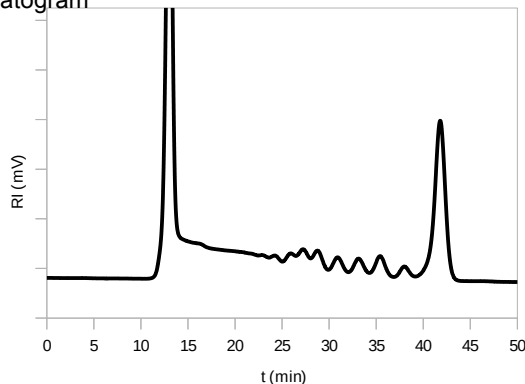



Detail enlargement

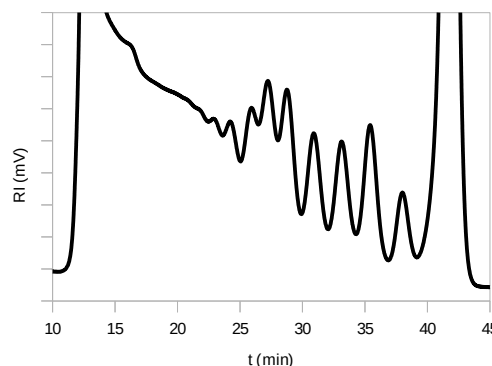

**AppliChrom ABOA SugarSep-Oligo-Ag, 300x8mm, Order No. SASOLAg103008**  
 "Maltodextrin 12 without extra glucose", Eluent: H<sub>2</sub>O, 0,25ml/min, 80°C

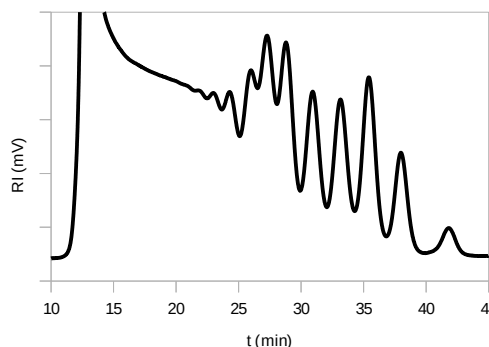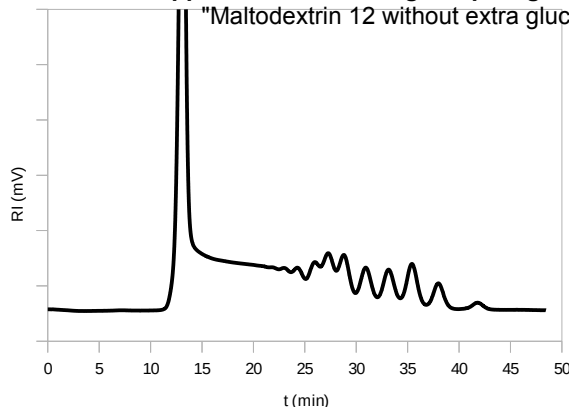

**For reason of Comparison: AppliChrom ABOA SugarSep-Oligo-Na, 300x8mm, Order No. SASOLNa103008**  
 "Maltodextrin 6", Eluent: H<sub>2</sub>O, 0,25ml/min, 80°C

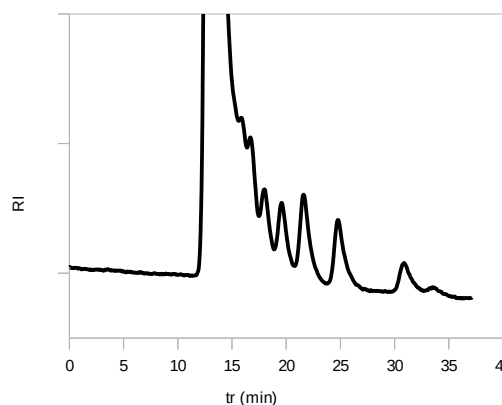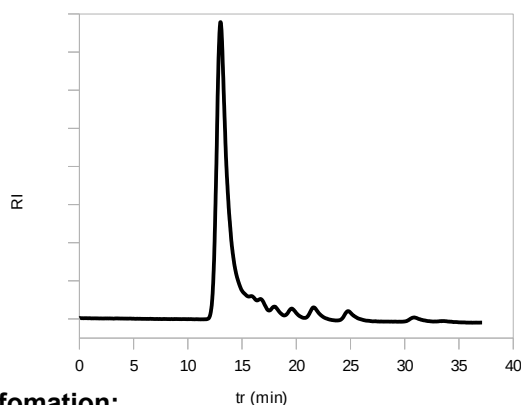

### Order Information:

SASOLAg101508	HPLC-Column AppliChrom® ABOA SugarSep-OligoAg, 10µ 150mm x 8mm	Price available upon request
SASOLAg103008	HPLC-Column AppliChrom® ABOA SugarSep-OligoAg, 10µ 300mm x 8mm	Price available upon request
SASOLAg10308	Precolumn AppliChrom® ABOA SugarSep-OligoAg, 10µ, 30mm x 8mm	Price available upon request
SASOLNa101508	HPLC-Column AppliChrom® ABOA SugarSep-OligoNa, 10µ 150mm x 8mm	Price available upon request
SASOLNa103008	HPLC-Column AppliChrom® ABOA SugarSep-OligoNa, 10µ 300mm x 8mm	Price available upon request
SASOLNa10308	Precolumn AppliChrom® ABOA SugarSep-OligoNa, 10µ, 30mm x 8mm	Price available upon request